

SystemClean mti-Hygiene Spray

Alcoholic rapid disinfection



Product description

mti-Hygiene Spray is a ready to use spray for alcoholic rapid disinfection of surfaces without reffatin agents. This product has a broad activity spectrum.

mti-Hygiene Spray is characterized by high disinfection performance on all alcohol-resistant, washable and wet-treatable objects and surfaces.

Wirkung

mti-Hygiene Spray is a special rapid and reliable surface disinfection eliminates 99.9% of bacteria, fungi, germs and certain viruses.

Our plus:

- Rapid
- Aldehyde free
- high disinfection performance

Application note

- clear the surface of visible contamination.
- Spray surfaces and objects until complete wetting with **SystemClean mti-Hygiene Spray**
- Leave on for at least 30 seconds. Leave on for 3 minutes for maximum effect. This process is repeated as required.

BAuA Registrier-Nr.: N88441

Active ingredients:

Ethanol, Isopropanol, fragrances -aldehyde free-

Store:

Room temperature

Material compatibility:

Suitable for alcohol-resistant surfaces.

Disposal:

Dispose of container to an approved disposal company.

Safety note:

Avoid contact with skin and eyes.

Order information

Cat. no.: 601-002 (1 x 1 L) Spray Bottle

Cat. no.: 601-002 (1 x 250 ml) Spray bottle with finger sprayer

